

Product use and handling instructions

Avoid contact among the magnets. In the case that two or more magnets contacted each other, do not pull the magnets apart; twist the magnets gently to prevent damage.

DRP-MagneTubes can be cleaned by spraying with 70% isopropyl alcohol and wiped with a dry cloth/paper. Avoid prolonged exposure to water or aqueous solutions. **Warning:** Do not autoclave DRP-MagneTubes.

Take special care when handling these magnets in the presence of magnetic devices that can be damaged by the magnetic field. In case of pacemaker users (or similar medical devices), the handling of this product is strongly not recommended.